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## GRANULATOR/ COMMINUTION EQUIPMENT

# NS

赛德力离心机制造专家  
SAIDELI CENTRIFUGE MANUFACTURING EXPERTS

www.saideli.com

**Saideli赛德力**

江苏赛德力制药机械有限公司  
JIANGSU SAIDELI PHARMACEUTICAL  
MACHINERY CO.,LTD.

制粒/粉碎设备  
GRANULATOR/COMMINUTION EQUIPMENT

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#### ● Introduction

YK-160A rock granulator is the equipment of prill and disintegrating with bottom discharging.

Structure and principle: The decelerating equipment is directly connected to the motor.

The eccentric cam power the rack, make the scraping axle spinning, and the solid-state materials into the inner-chamber of the scraping axle, then the materials are push out of the screen mesh of the point angle of the axle in the shape of solid prills or powder. Distinguishing feature: This granulator is very capable for prill, Its structure is simple, the running is well, and the repairing is convenient. Its volume is small.

Main usage: it is mainly used in the trade of pharmacy and chemical foodstuff. And is suitable for prilling the wet powder, smashing the dried power.

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#### ● Introduction

This machine is the necessary equipment in the trade of chemical and pharmacy for the shiver of solid state materials the emulsification of liquid materials.

Work principle: The movement, such as impact, collision to the materials from the special high-speed rotor cause the high-speed movement of the materials inside the machine. The course of the movement is the course of the sieving and smashing.

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